IMMB-09-02-05-22-EN-PL

Start up Guide

Szybki start

Precision Balances

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- PM C32

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
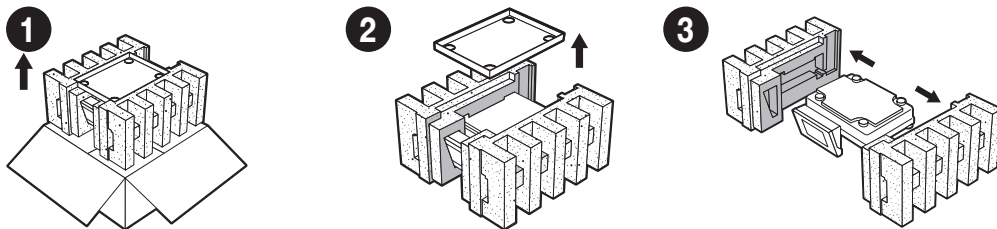
PM C32 (max: 10-50 kg)

PM C32 (max \geq 60 kg)

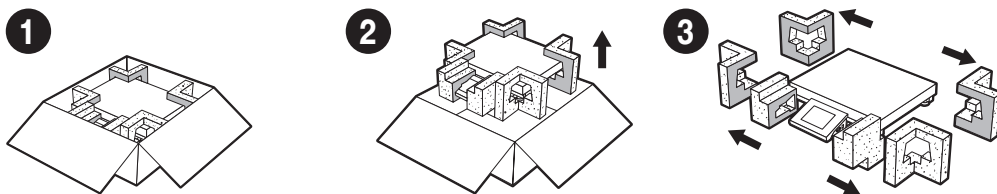
Keep the packaging should a warranty claim or service be required.

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3. RECOMMENDATIONS AND WARNINGS

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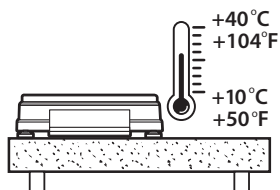
Operate the device in a room where the temperature ranges between 10–40 °C (50–104 °F) and where the relative humidity is below 80%. Temperature must change slowly and gradually.


Using the anti-vibration table is recommended in order to enable the most stable and repeatable measurement.

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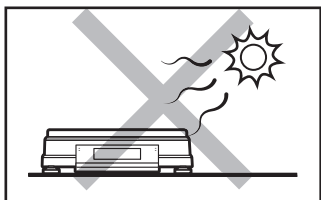
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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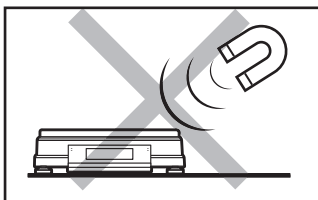
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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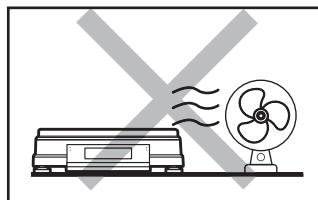
Avoid air drafts and air movements at the workstation.

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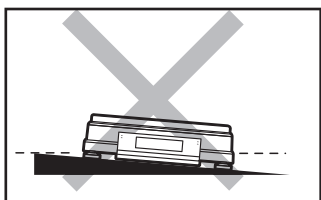
Make sure that the balance is placed on an even surface.

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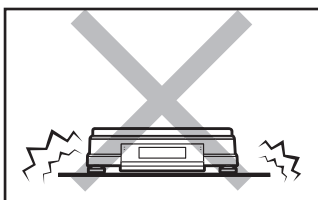Do not place the balance on an unstable surface exposed to shocks and vibrations.

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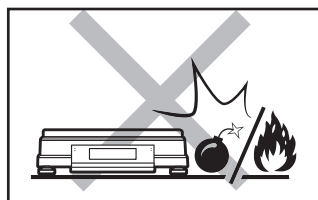Not intended for usage within EX zones. Not intended for weighing explosive or flammable substances.

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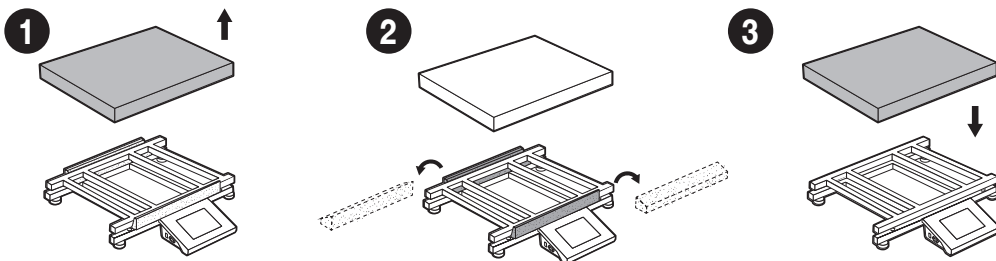
4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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4.1. Remove transport lock

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
PM C32 (max ≥ 60 kg)

4.2. Grounding spring check

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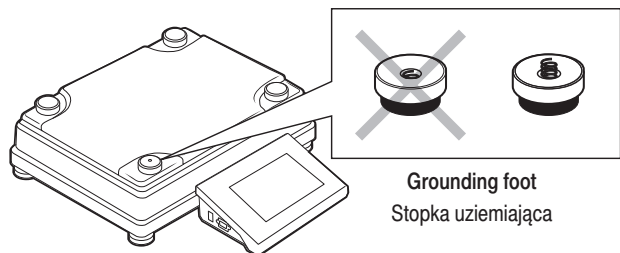PM C32 (max: 10-50 kg)


Check whether the grounding spring takes its intended location. Make sure that the grounding spring juts slightly out of the hole.

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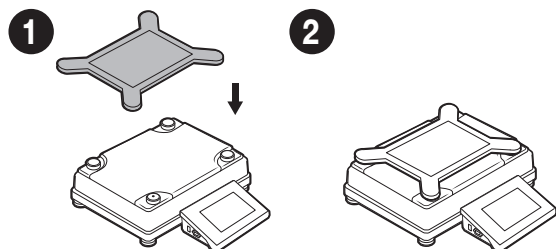
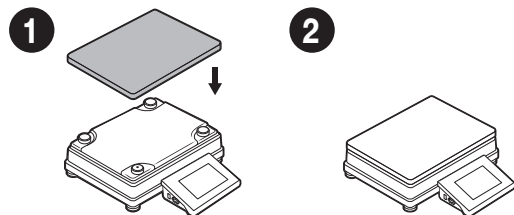
PM C32 (max ≥ 60 kg)

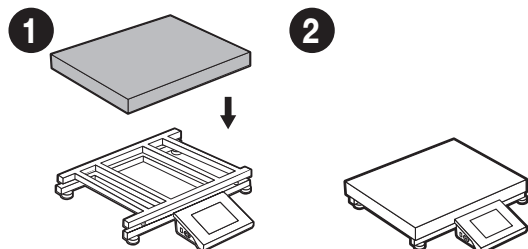6. PREPARING FOR WORK

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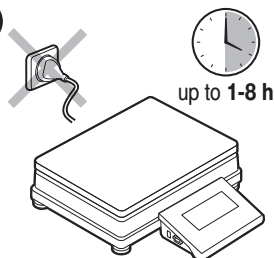
Place the device at the workstation, wait until it reaches room temperature.

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6.1. Temperature stabilization

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
1

Prior to switching the device on, let it reach room temperature (estimated stabilisation time: 1-8 hours).

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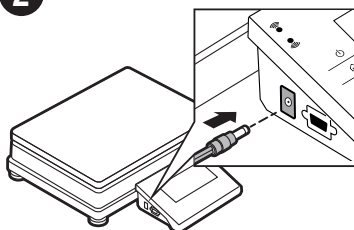
2

Connect the power adapter to DC connector.

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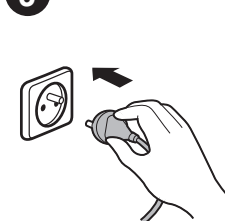3


Connect the power adapter to the mains.

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6.2. Balance levelling

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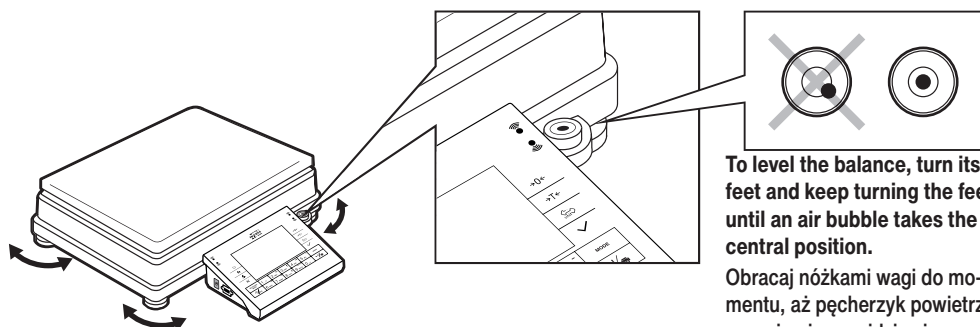
PM C32 (max: 10 - 50 kg)

To level the balance, turn its feet and keep turning the feet until an air bubble takes the central position.

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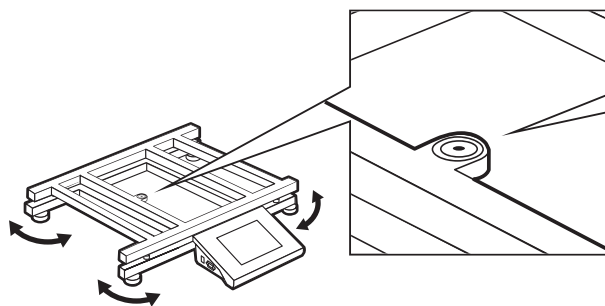
6.3. Switching the balance on

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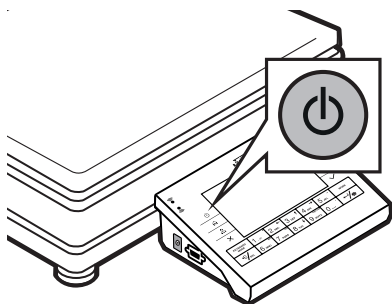To switch the device on, press ON / OFF key.

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

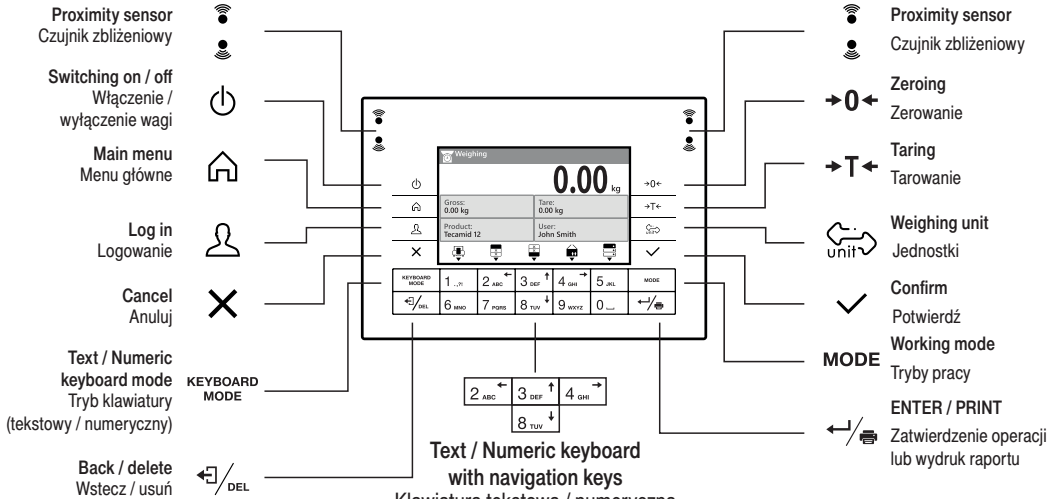
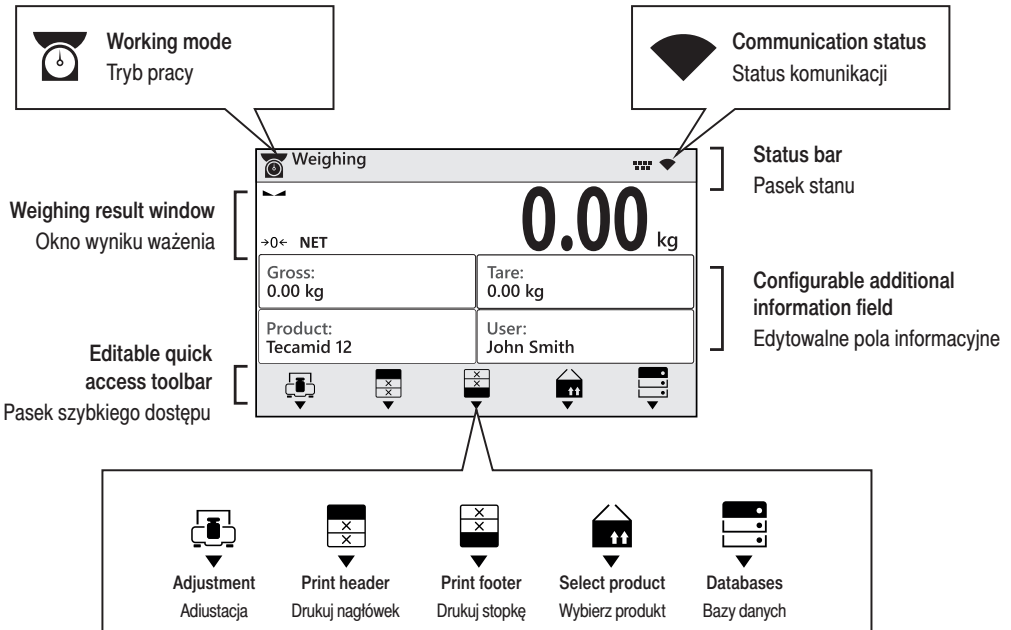
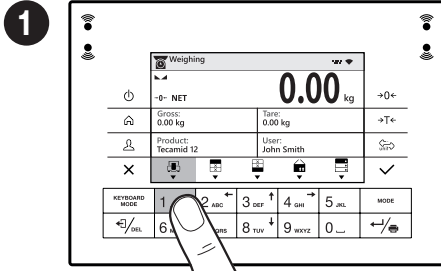
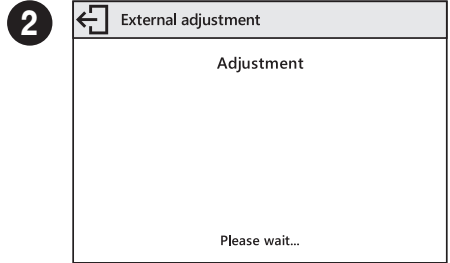
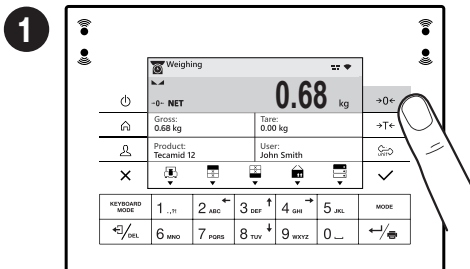
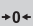
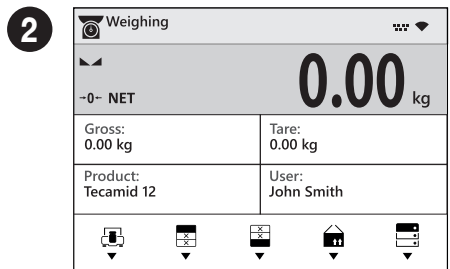
8.3. Taring

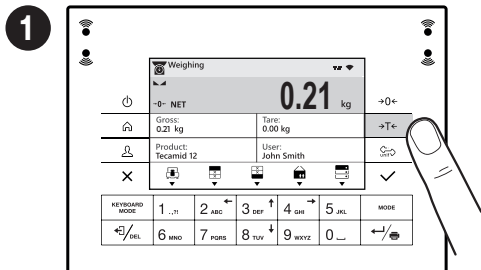Tarowanie

With loaded weighing pan: upon stabilization of the weighing result, press **→T←** key.

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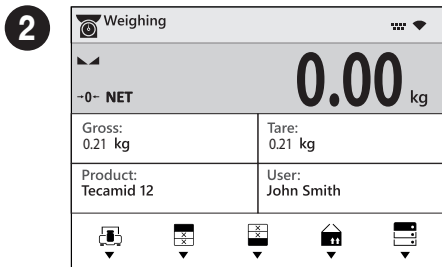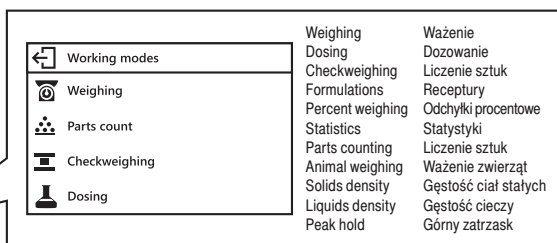
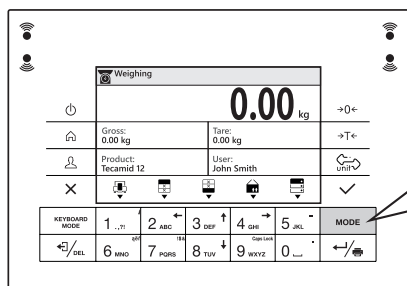
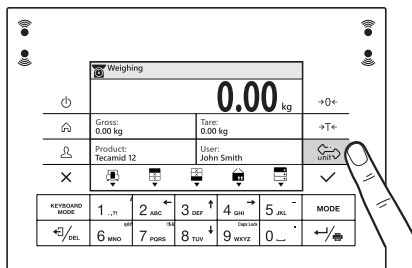
8.5. Units selection

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Press the **↵** key to change the weight unit.

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The balance enables selecting the following weighing units:

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9. SETTINGS

USTAWIENIA

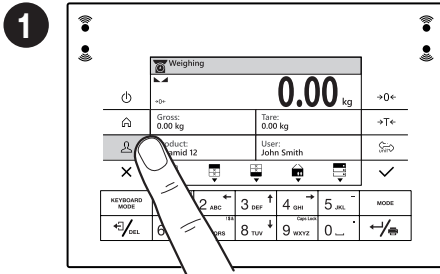
Some balance settings are accessible for Administrator exclusively.
Prior balance parameters setup, log in as the Administrator.

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

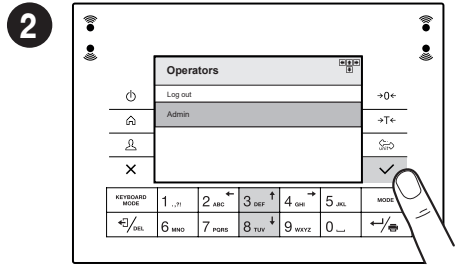
Users list is displayed. Using   keys, select "Admin" option and press  key to confirm.

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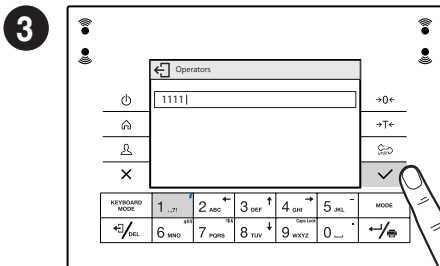Using numeric keyboard enter "1111" password and press  key to confirm.

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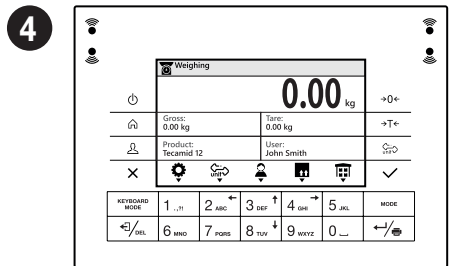9.2. User settings

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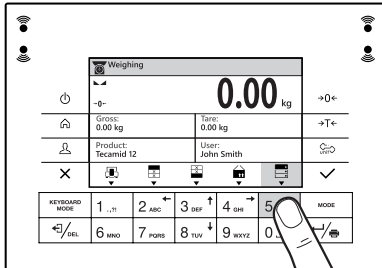
Prior balance users setup, log in as Administrator (point 9.1).

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
1

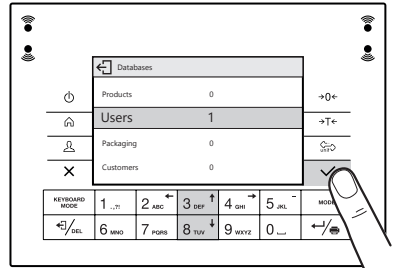Press the key under the  icon to enter databases.

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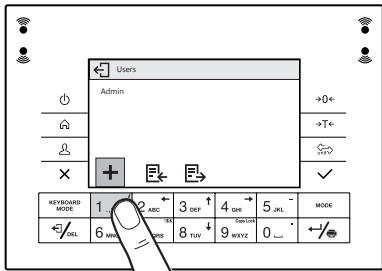
2

Using   keys, go to “Users” and press  key to confirm.

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
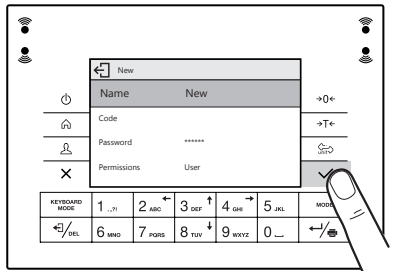
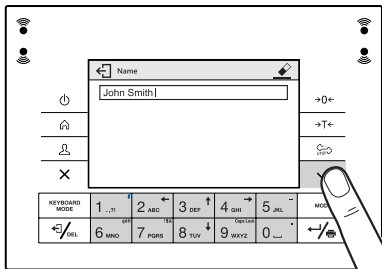
5

Using keyboard enter user name and press  key to confirm.

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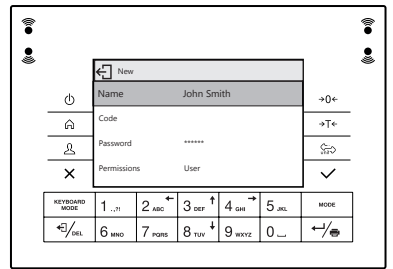7 Following the above procedure, set remaining parameters:

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Password: Enter user password

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Access levels: Set respective access level (user, advanced, admin)

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New	
Name	John Smith
Code	
Password	*****
Permissions	User

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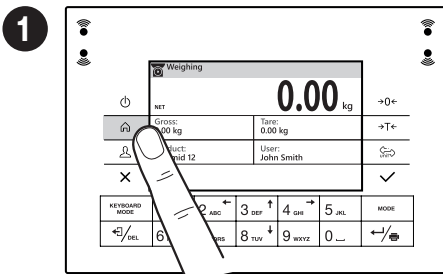
Prior balance proximity sensors setup, log in as Administrator (point 9.1).

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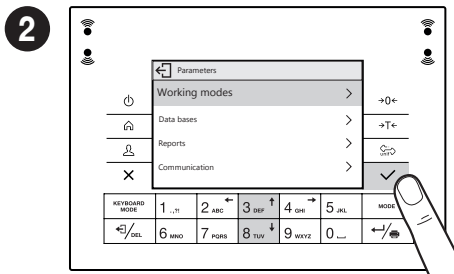
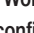
You can assign the proximity sensors with different functions for each working mode (the sensors will trigger different operation for each working mode). The diagram presents sensors setup for the “Weighing” mode.

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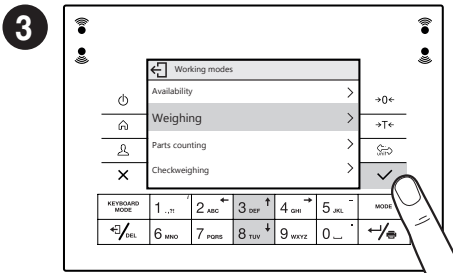
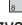
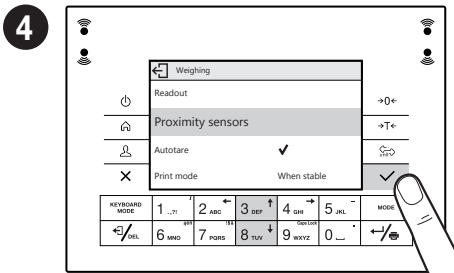
Using  keys, go to “Working Modes” and press  key to confirm.

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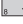
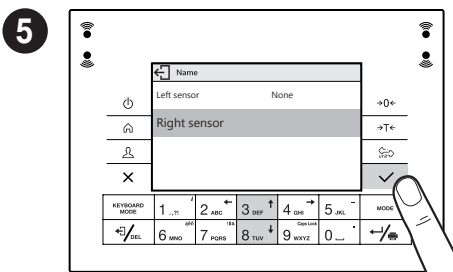
Using   keys select “Weighing” mode, and press  key to confirm.

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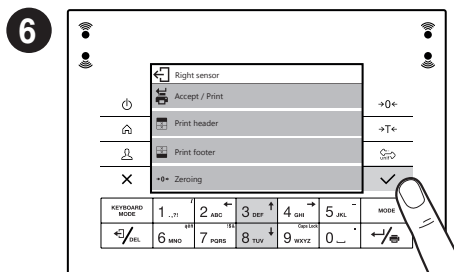
Using   keys select “Proximity sensors”, and press  key to confirm.

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Using   keys select the appropriate sensor (right or left), and press  key to confirm.

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Using   keys set function that is to be assigned to a selected sensor, and press  key to confirm.

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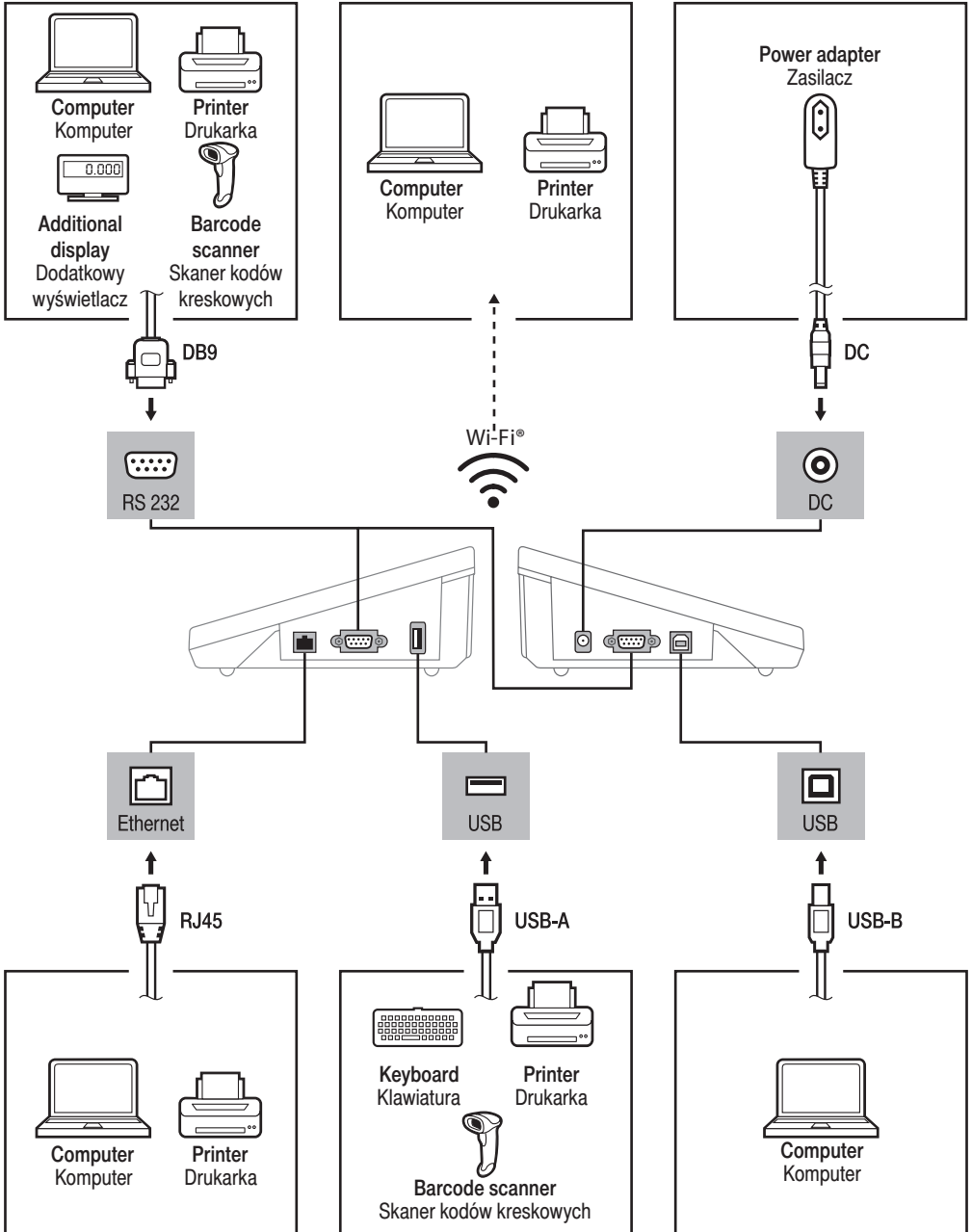
Following the above procedure you can configure sensors for other working modes.

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Wi-Fi® is a registered trademark of Wi-Fi® Alliance.

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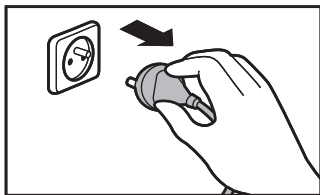
11. DEVICE CLEANING

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
11.1. General recommendations

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
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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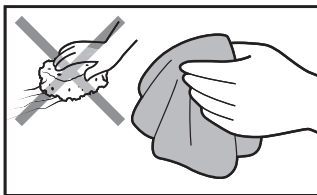Use soft cloths made of microfiber, natural fiber or man-made fiber. Avoid using abrasive cloths or cloths that might scratch the surface.

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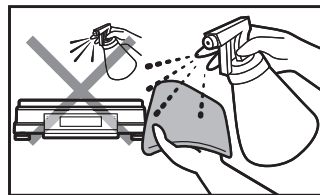Apply the cleanser onto the cloth first. Avoid applying the cleanser directly onto the device.

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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