IMMB-14-02-03-22-EN-PL

Start Up Guide

Szybki start

Ultra-Microbalances and Microbalances

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
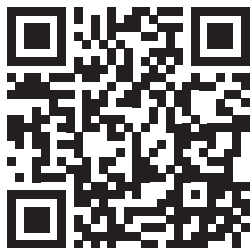
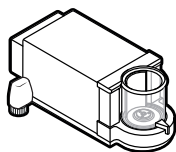
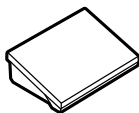
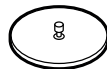
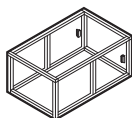
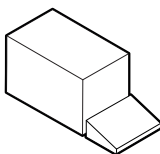
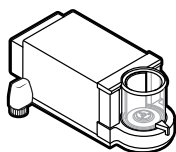
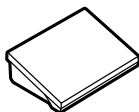
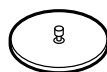
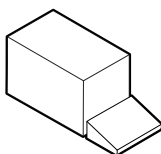
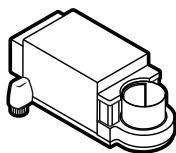
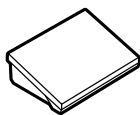
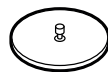
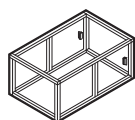
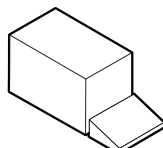
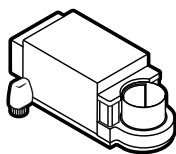
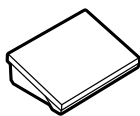
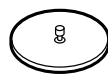
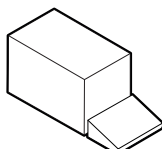
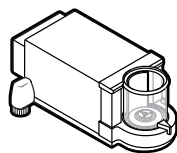
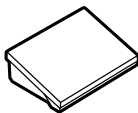
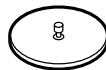
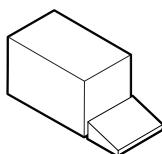
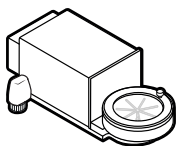
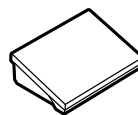
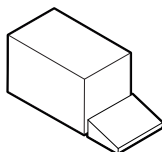
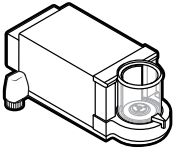
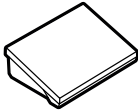
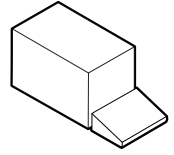
x1

2. UNPACKING

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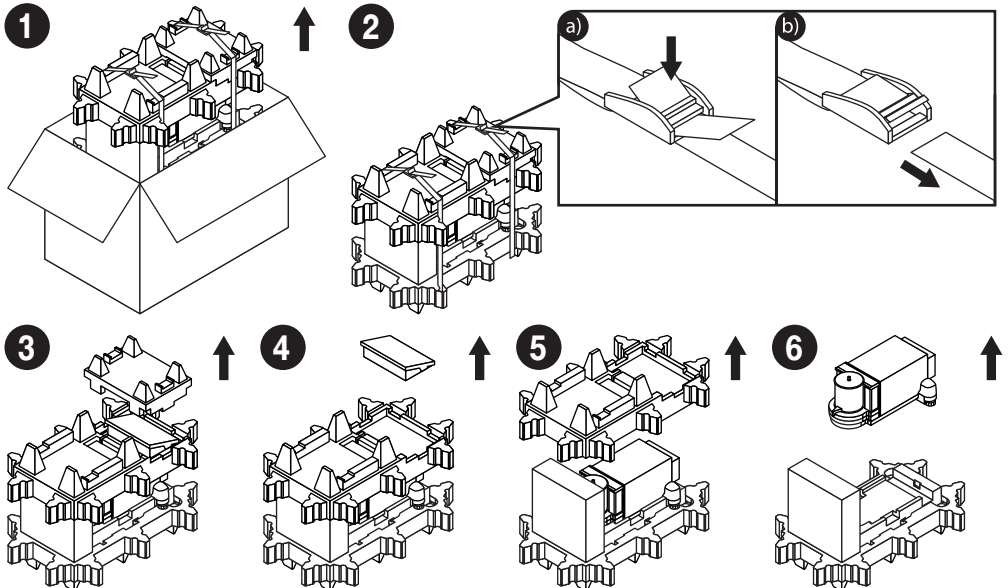
Keep the packaging for warranty and service purposes.

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3. RECOMMENDATIONS AND WARNINGS

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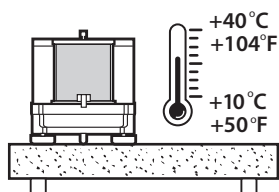
Operate the device in a room where the temperature ranges between 10 – 40 °C (50 – 104 °F) and where the relative humidity is below 80 %. Place the balance on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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
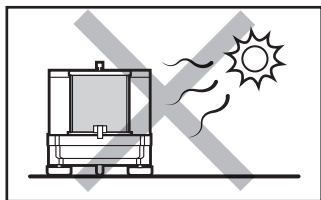
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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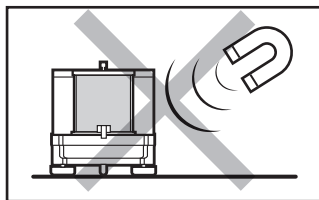
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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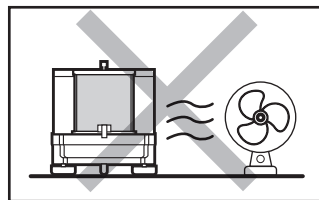
Avoid air drafts and air movements at the workstation.

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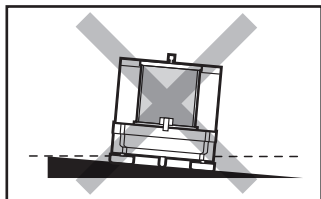
Make sure that the balance is placed on an even surface.

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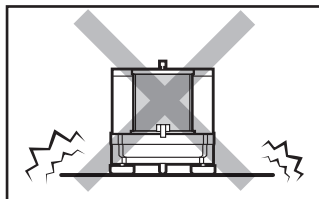Do not place the balance on an unstable ground exposed to shocks and vibrations.

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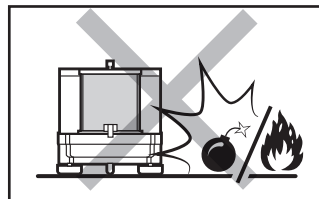
Do not place the balance in a hazardous area. Do not weigh explosive materials.

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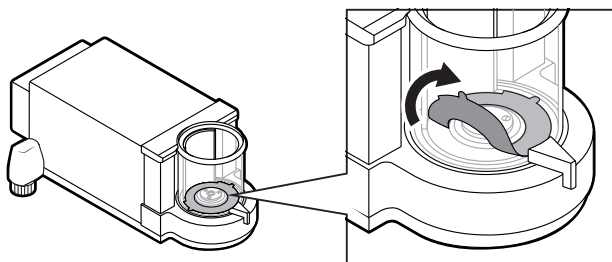
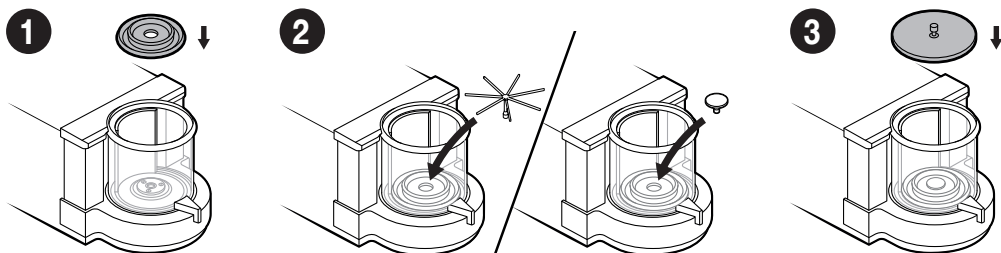
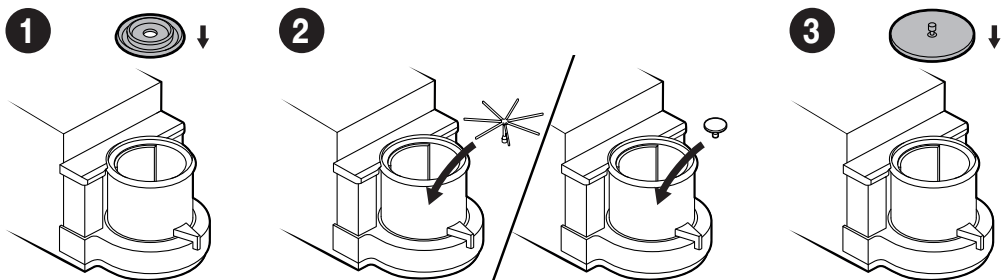
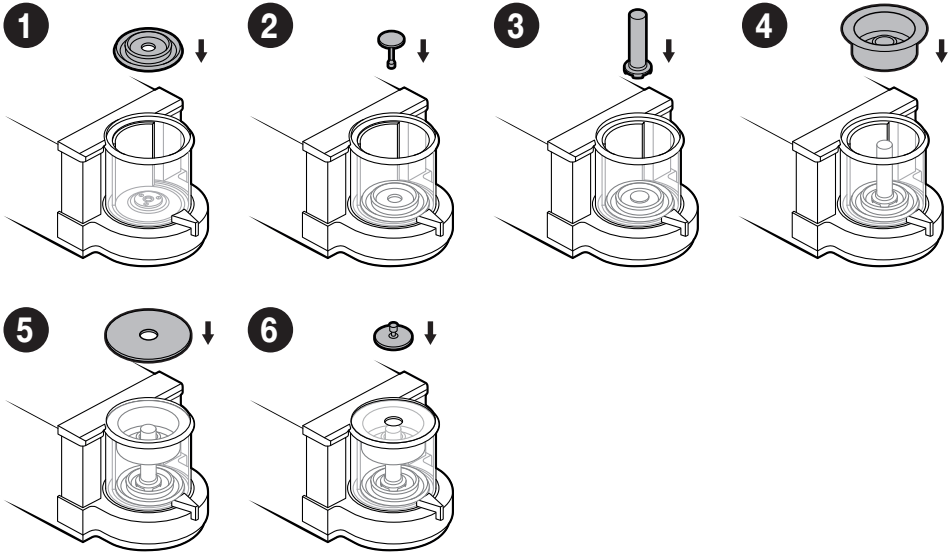
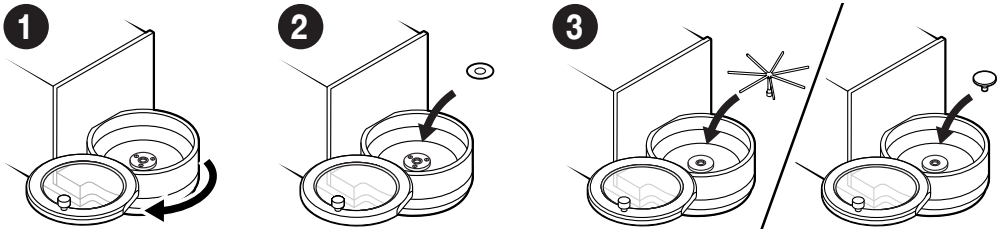
5.2. Connecting the Terminal

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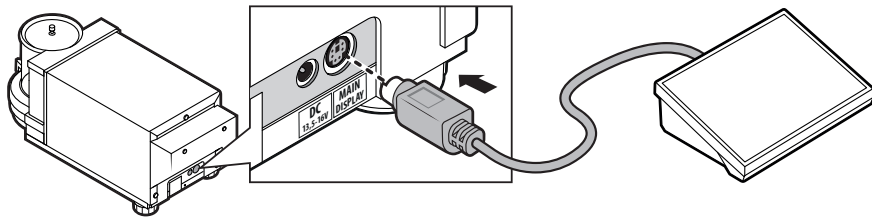6. PREPARING FOR WORK

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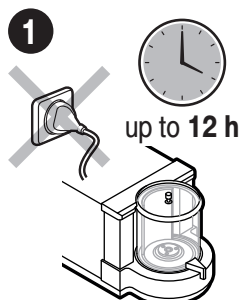
Place the device at the workstation, wait until it reaches room temperature.

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
6.1. Temperature Stabilization

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
Prior to switching the device on, let it reach room temperature (estimated stabilization time: up to 12 hours).

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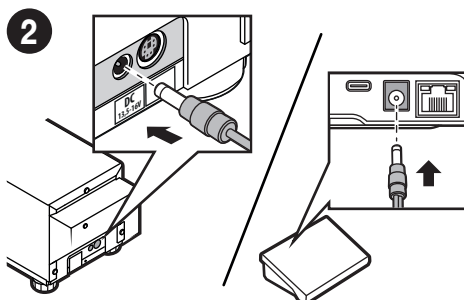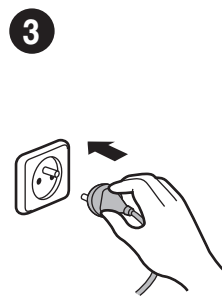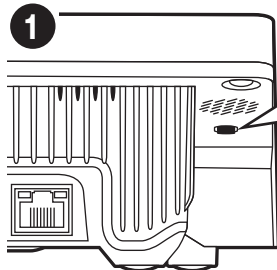
Do not connect the balance and the terminal to the power supply at the same time.

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
6.2. Switching the Balance On

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The balance switches on automatically when connected to the mains. If the balance is manually switched off using the top menu, it must be switched on using the button underneath the display housing on the left.

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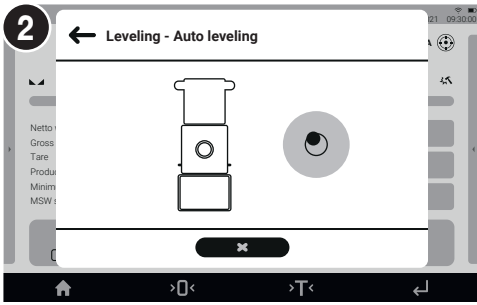
6.3 Manual Levelling

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
Ultra-microbalances and microbalances are equipped with the Reflex Level System, which monitors and automatically corrects the balance level.

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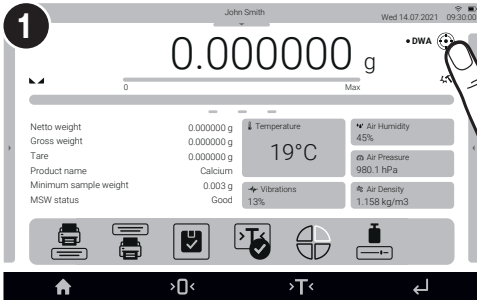
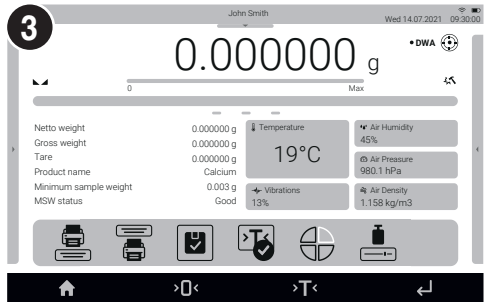
When autolevelling is complete, the home screen is displayed.

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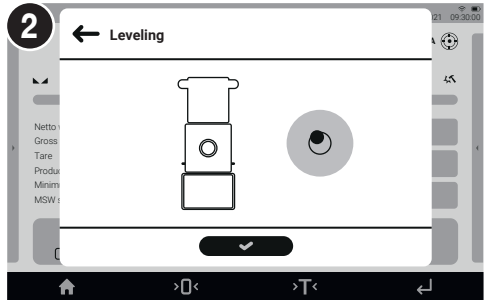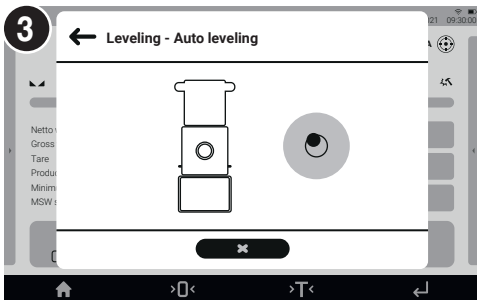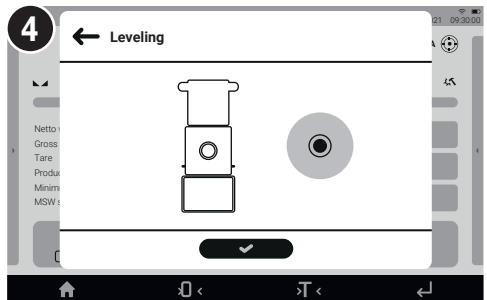
7. HOME SCREEN

EKRAN GŁÓWNY

information about the currently used working mode, logged-in operator, date, time, active connection with a computer.

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The screenshot shows the Home Screen with the following data:

- Netto weight: 0.000000 g
- Gross weight: 0.000000 g
- Tare: 0.000000 g
- Product name: Calcium
- Minimum sample weight: 0.003 g
- MSW status: Good
- Temperature: 19°C
- Air Humidity: 45%
- Air Pressure: 980.1 hPa
- Vibrations: 13%
- Air Density: 1.158 kg/m³

Additional information shown includes: Weighing Fast Dosing, John Smith, Wed 14.07.2021 09:30:00, and a DWA status indicator.

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8. BASIC OPERATIONS

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8.1. Internal Adjustment

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The screenshot shows the Home Screen with the Internal Adjustment button (represented by a house icon) highlighted with a red circle and the number 1. The rest of the screen content is identical to the previous screenshot.

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
The screenshot shows the Adjustment menu with the following options:

- Adjustment
- Operators
- Profiles
- Database
- Communication
- Peripherals
- Administrator panel
- Miscellaneous
- Ambient conditions
- Scheduled tasks
- Update
- About

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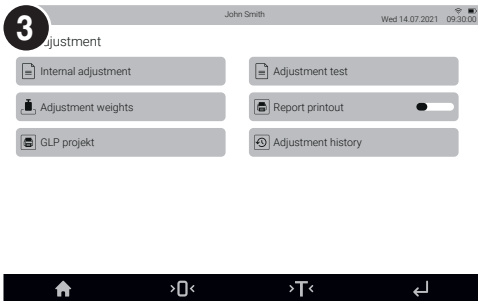
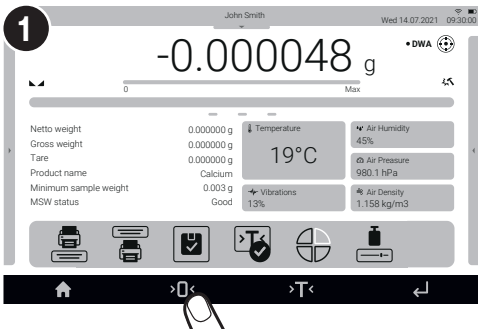
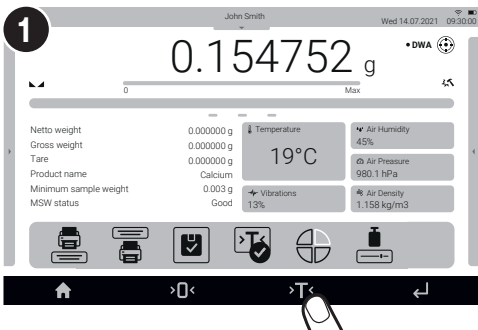
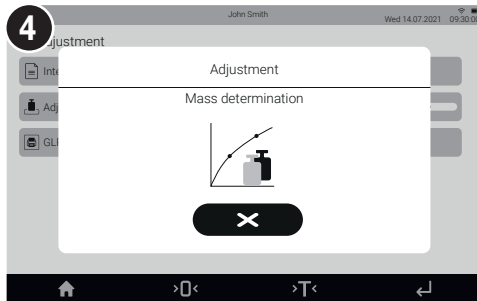
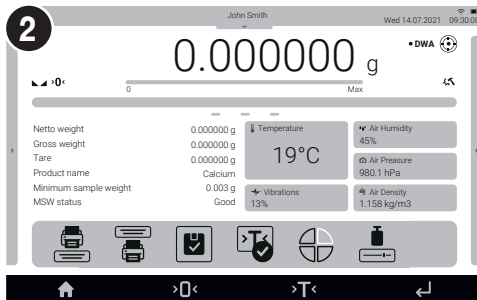
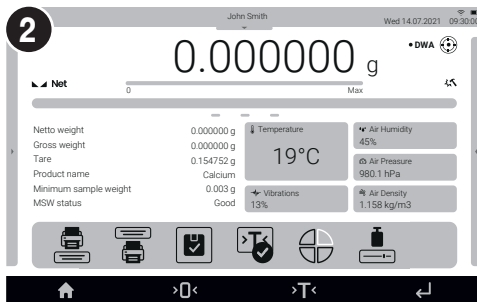
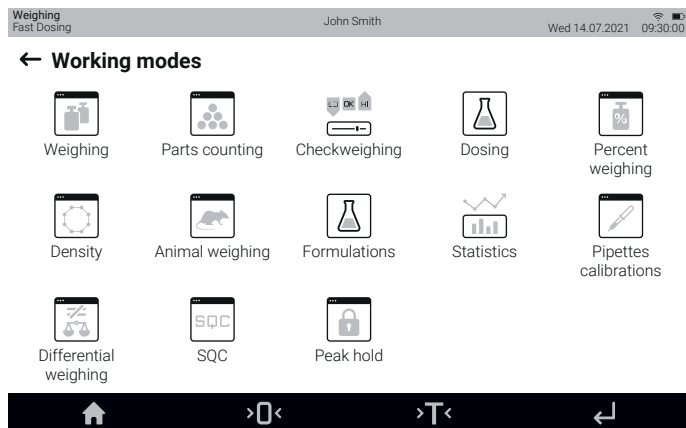
8.5. Unit Selection

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Press the current mass unit pictogram to display the list of available units.

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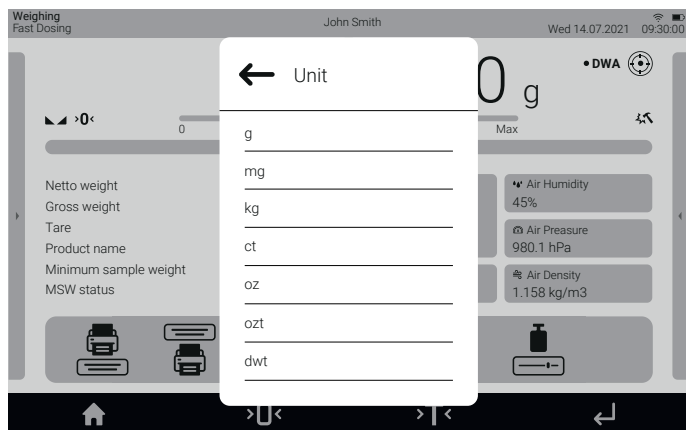9. SETTINGS

USTAWIENIA

Some balance settings are available only for the Administrator.
Prior to the balance parameters setup, log in as the Administrator.

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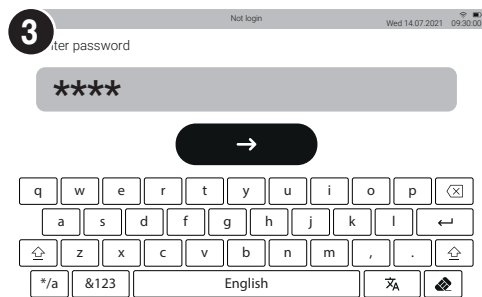
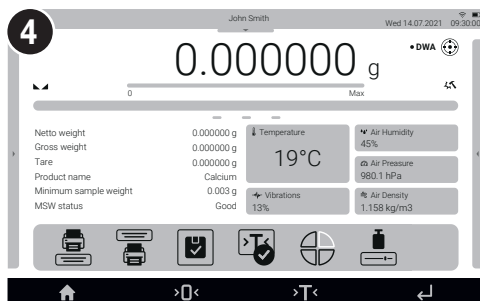
You have logged as the Administrator.

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9.2. Operator Settings

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Prior to operators configuration, log in as the Administrator (see section 9.1).

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
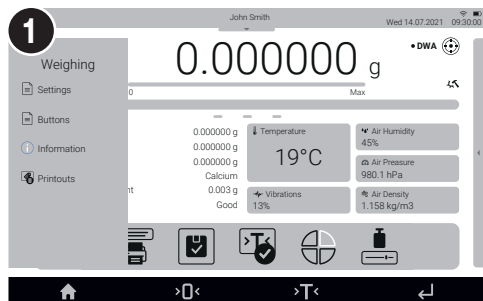
You can assign a function of your choice to each proximity sensor. In addition, you can assign various functions to the same sensor in different working modes. See the exemplary procedure for sensors configuration for the “Weighing” mode.

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Slide out the menu.

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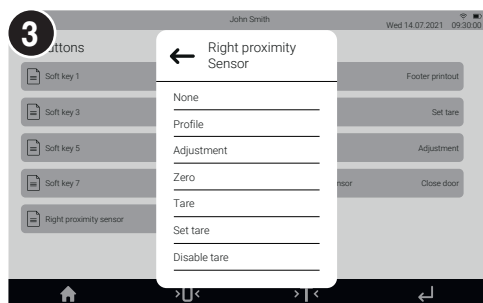
Select a respective sensor (right or left) from the list.

Select a function to be assigned to the sensor.

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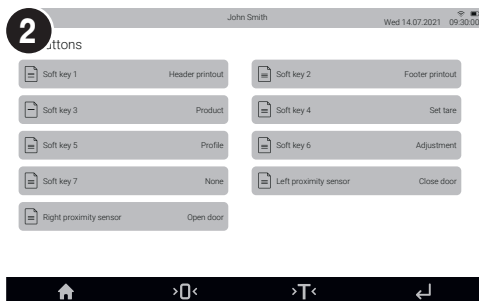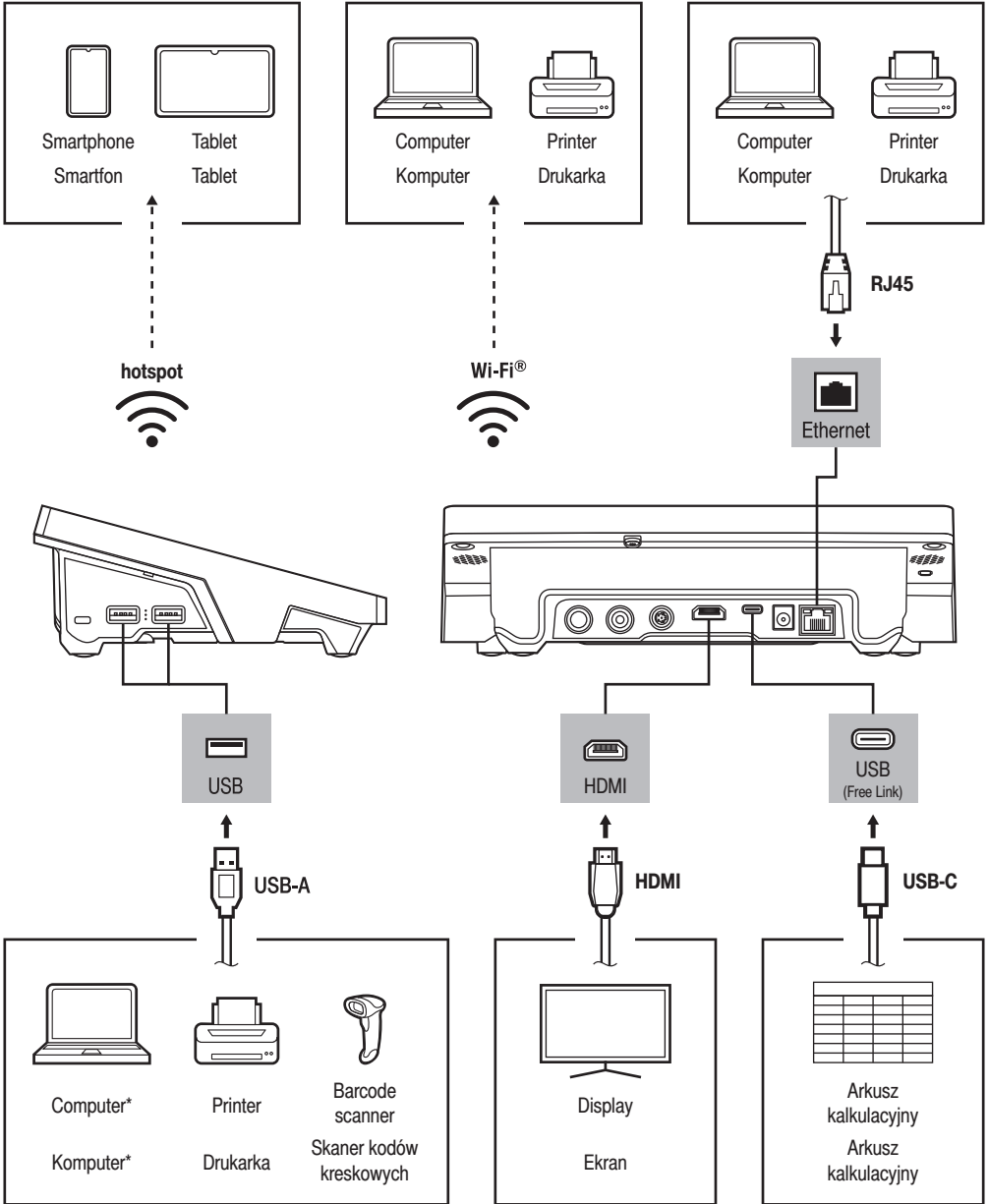
Wi-Fi ® is a registered trademark of Wi-Fi Alliance.

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*A USB to RS232 converter must be used for proper connection.

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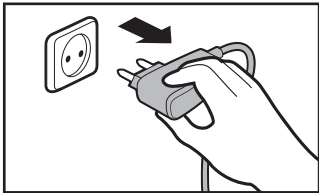
11. CLEANING AND MAINTENANCE

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
11.1. Guidelines and Precautions

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
Prior to cleaning, it is necessary to disconnect the device from the mains! Remove the plug from the outlet and disconnect all cables.

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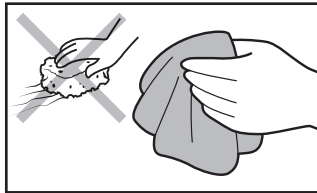Use soft cloth made of either natural or synthetic fibers exclusively. Avoid using stiff or hard fabric, as this may cause scratches on the surface!

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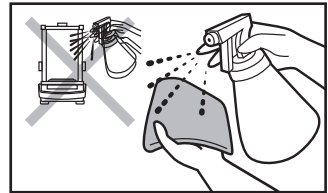Apply the cleanser to the cloth. Remember not to apply the cleanser directly onto the device!

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Use neither sharp-structured nor scrubbing-purposed substances.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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If it is necessary to disassemble the weighing chamber, follow the instructions in section 11.3.

- **GLASS PANES AND COMPONENTS** can be cleaned using window cleaner.
- **WEIGHING PANS AND STAINLESS STEEL OR ALUMINIUM COMPONENTS** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- **OPERATION PANEL AND HOUSING** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- Dry sample leftovers can be removed using either a brush or a handheld vacuum cleaner.
- Clean components can be dried using a soft cloth or a dust-free paper towel. With this, the remaining moisture will be absorbed.
- **Install all the components, providing they are completely dry.**

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
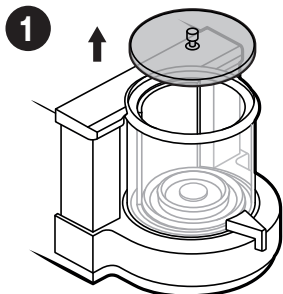
Prior to disassembling the weighing chamber elements, make sure the microbalance is turned off and unplugged from the mains. When disassembling the elements, use the tweezers that come standard with the instrument.

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
Lift the lid of the weighing chamber.

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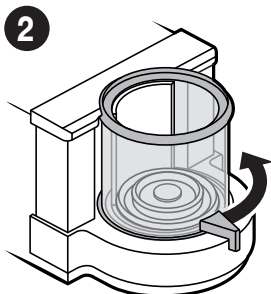
Open the weighing chamber by twisting the handle.

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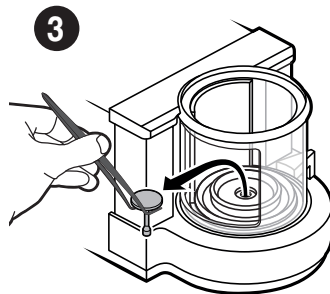
Using the tweezers, carefully remove the weighing pan from the weighing chamber.

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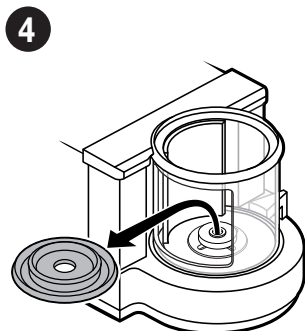
Remove the draft shield from the weighing chamber.

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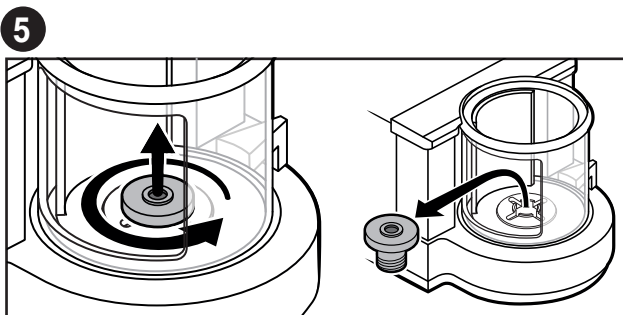
Unscrew the draft shield's sleeve counterclockwise and remove it from the weighing chamber.

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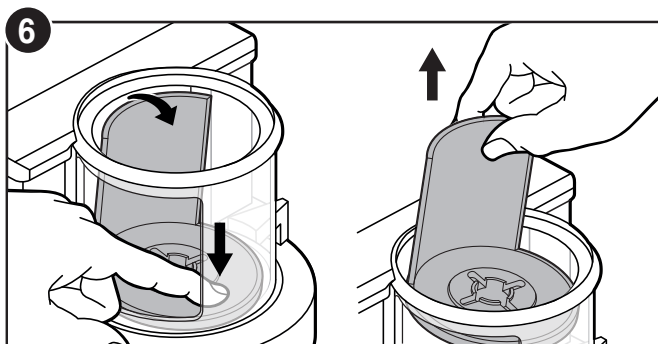
6 Use your finger to press the base of the ring assembly in order to tilt the top edge. Grab the top edge and remove the ring assembly through the upper hole of the weighing chamber.

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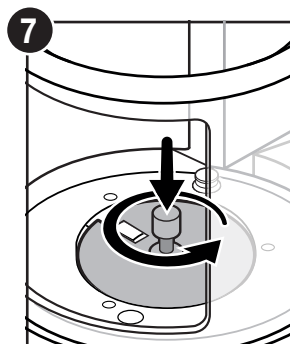
7 Install and tighten the protective ring using the handle.

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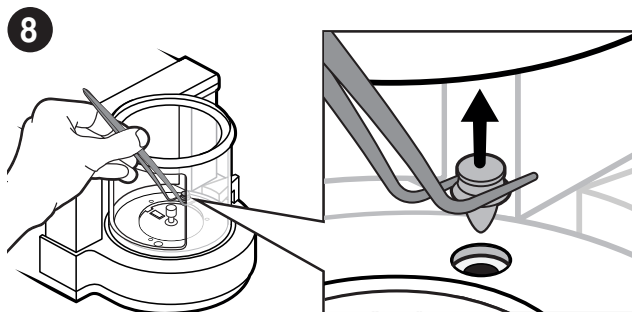
8 Remove the plastic fastener from the weighing chamber using tweezers.

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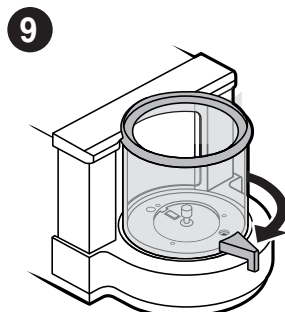9 Close the weighing chamber.

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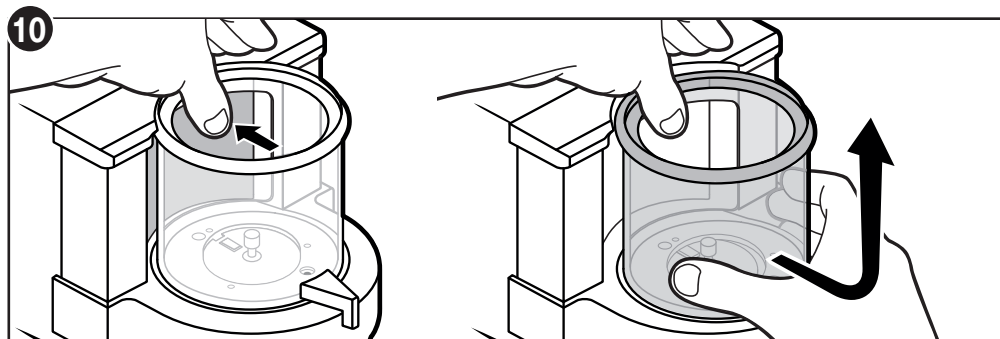10 Press the weighing chamber clamp. While pressing the clamp, grab the chamber, move it towards you and up.

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After cleaning the weighing chamber elements, they must be assembled in the reverse order.

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