ITKB-10-02-05-22-EN-PL

Start up Guide

Szybki start


Indicator

Miernik

PUE 7.1

To view the full user manual, go to the website or scan the QR code:

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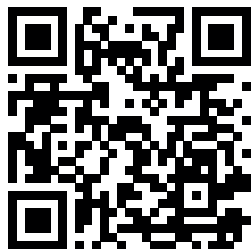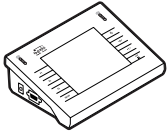
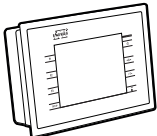
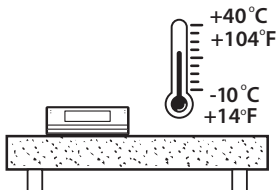
2. NOTES AND WARNINGS

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
Operate the indicator in a room free from dust where the temperature ranges -10 – 40 °C ($+14$ – 104 °F) and where the relative humidity is below 80%. Potential temperature change shall occur gradually and slowly.


Place the scale on a table or a wall bracket ensuring stability. For the most stable and repeatable results it is recommended to use an anti-vibration table.

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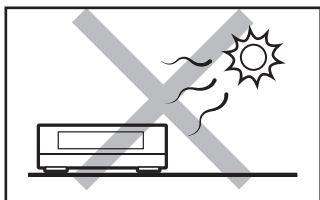
Place the scale away from heat sources. Avoid exposing the scale to the sunlight.

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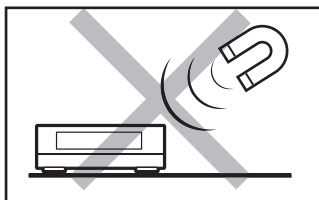Avoid exposing the scale to the influence of a magnetic field. Do not weigh magnetic substances.

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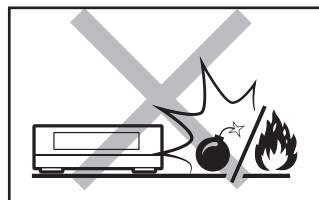Not intended for use within EX zones. Not intended for weighing explosive or flammable substances.

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3. PREPARING FOR WORK

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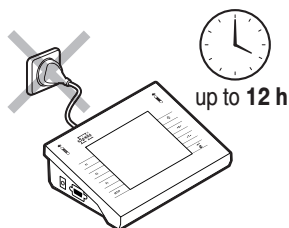
Place the indicator on the workstation and leave it until it reaches ambient temperature.

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3.1. Indicator Temperature Stabilization

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
1

Prior to switching the indicator on let it reach room temperature (estimated stabilization time: 12 hours).

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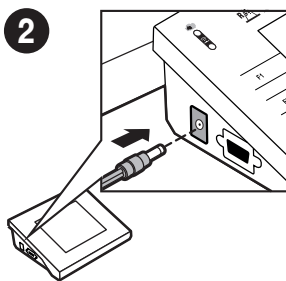
2

Connect the power supply to a connector located on the side of the indicator.

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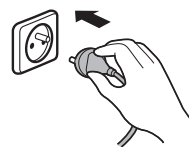3


Connect the power supply to the mains.

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3.2. Indicator Start-Up

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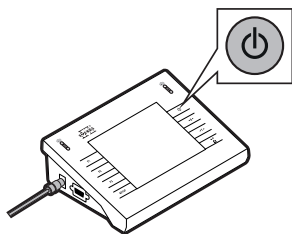Press ON / OFF key to start the indicator.

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The indicator is ready for operation!

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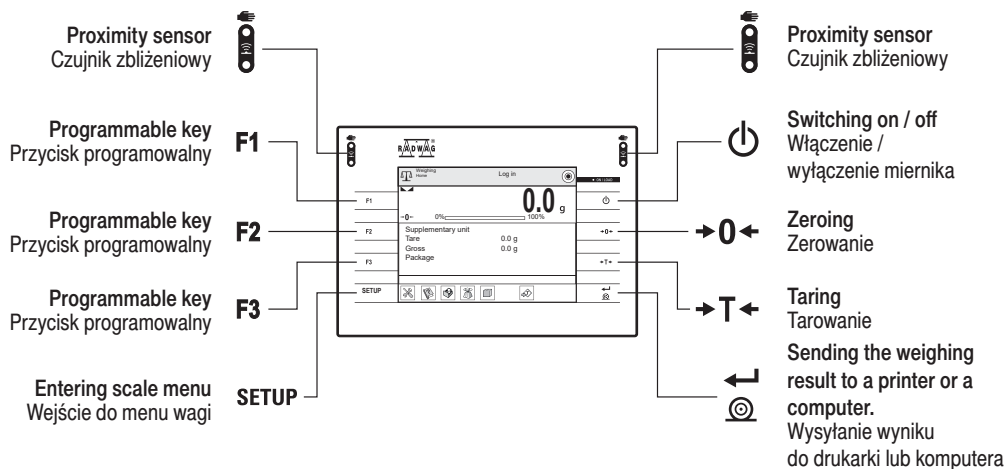
5.1. Working Modes / Unit Selection

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Working modes

1	Weighing
2	Parts counting
3	Percent weighing
4	Dosing
5	Formulations
6	PGC

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

5.2. Working Modes Settings

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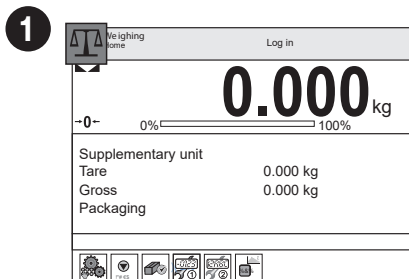
Press the current mass unit pictogram to display the list of all available units.

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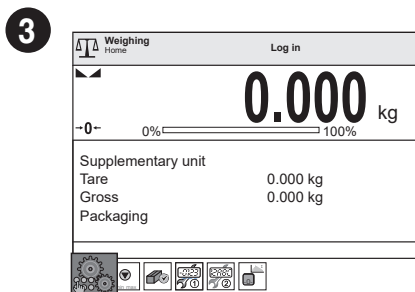
Press the working mode settings pictogram to edit the current working mode.

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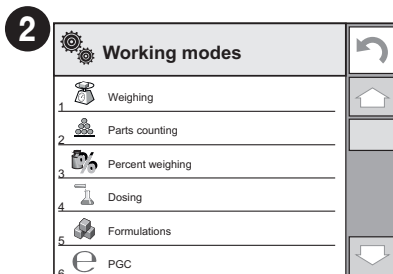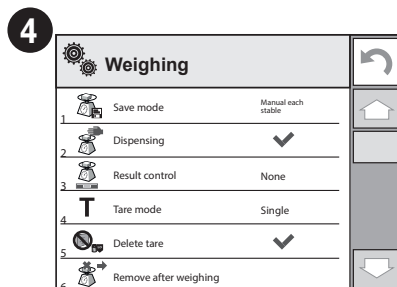A list of editable parameters of a given working mode is displayed.

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6. SETTINGS

USTAWIENIA

Some scale settings are only available to the Administrator.

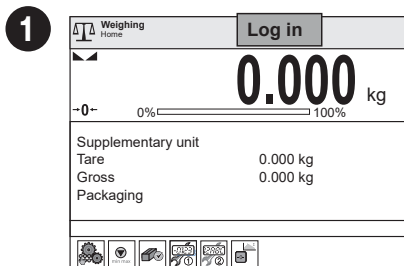
Prior to setting scale parameters log in as the Administrator.

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

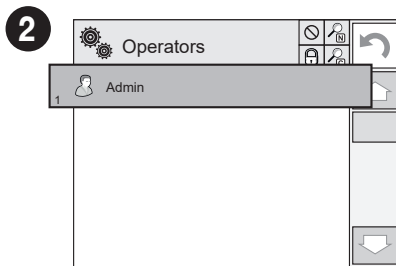
Select „Administrator” from the displayed list of operators.

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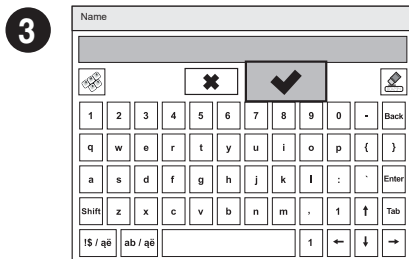When the field for filling the password is opened, confirm the empty field by pressing key.

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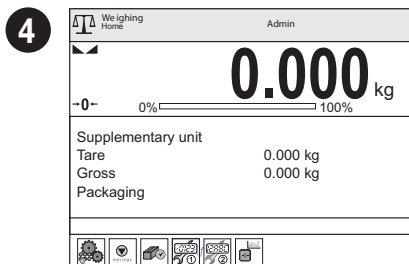You have logged as the administrator.

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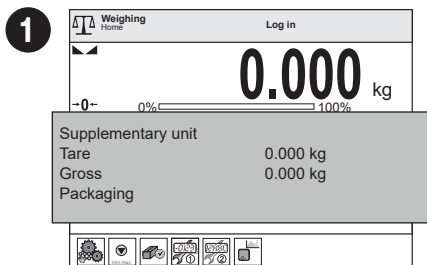
Prior to configuring the proximity sensors, log in as the Administrator (see step 6.1.).

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
You can assign a function of your choice to each proximity sensor. In addition, you can assign various functions to the sensor in different working modes. See the exemplary procedure for sensors configuration for the “Weighing” mode.

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
Press the workspace to access current working mode settings.

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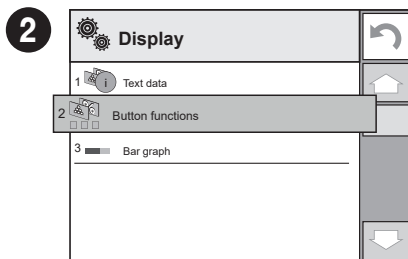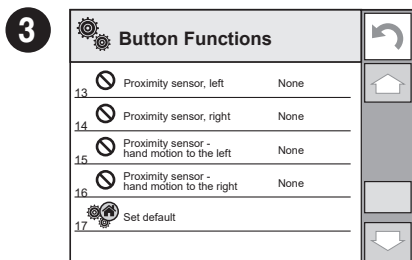
Select respective sensor (right or left) from the list.

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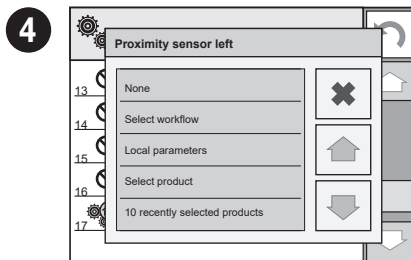Select function to be assigned to the sensor.

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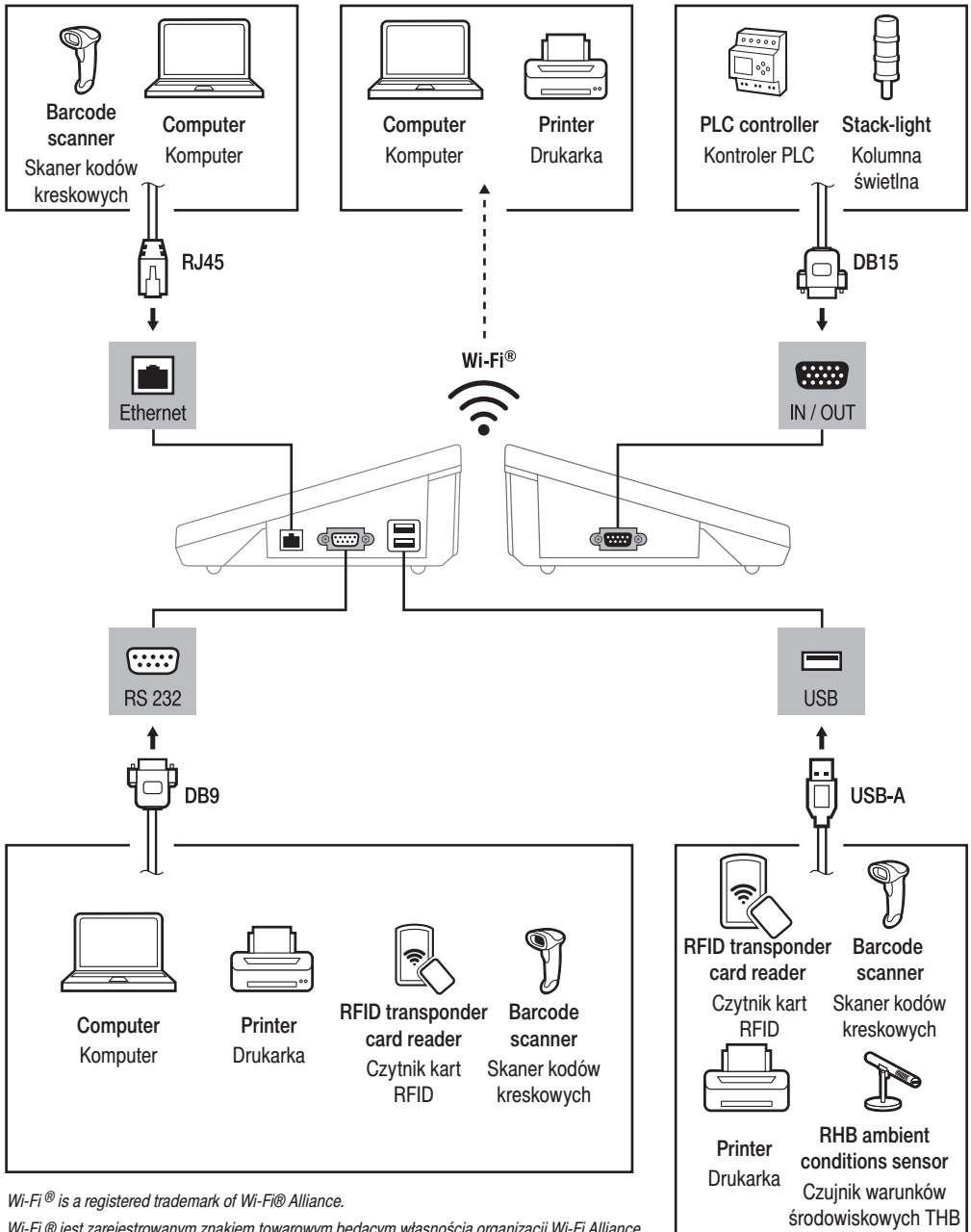
Follow the above procedure to configure proximity sensors for other working modes.

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Wi-Fi® is a registered trademark of Wi-Fi® Alliance.

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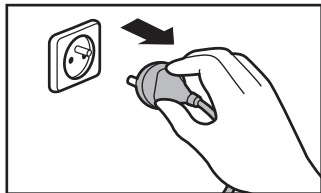
8. CLEANING AND MAINTENANCE

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8.1. Notes and Warnings

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
Prior to cleaning it is necessary to disconnect the indicator from the mains. Unplug the connector from the electrical socket.

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Avoid using aggressive cleaning agents (e.g. solvents, cleansers containing chlorine, corrosive substances and bleach).

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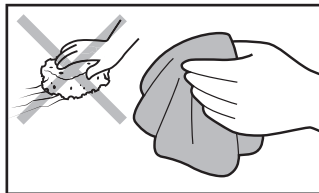
For cleaning use only soft cloths made of microfiber. Do not use any abrasive cloths or any material that may scratch the surface!

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Avoid using cleansers containing abrasive substances or scouring preparations.

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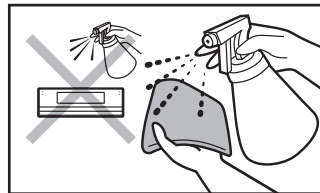Apply the cleanser onto the cloth first. Avoid applying the cleanser directly onto the device!

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Make sure that no dust or liquid enters the sockets and the inside of the indicator.

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