ITKB-08-02-05-22-EN-PL

Start up Guide

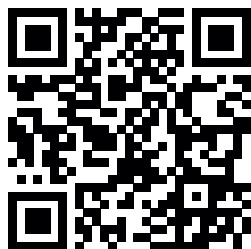
Szybki start

Indicator

Miernik

- PUE C315

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Indicator
Miernik
x1

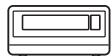
Power supply*
Zasilacz sieciowy*
x1

**The type of plug may vary depending on the country.*

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2. NOTES AND WARNINGS

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
+40 °C
+104 °F
-10 °C
+14 °F


Operate the indicator in a room free from dust where the temperature ranges -10–40 °C (+14–104 °F) and where the relative humidity is below 80%. Potential temperature change shall occur gradually and slowly. Place the scale on a table or a wall bracket ensuring stability. For the most stable and repeatable results it is recommended to use an anti-vibration table.

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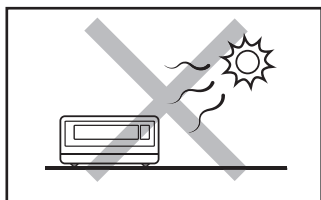
Place the scale away from heat sources. Avoid exposing the scale to the sunlight.

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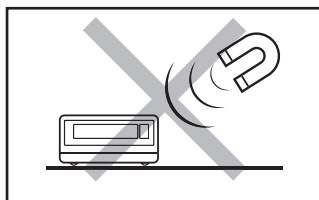Avoid exposing the scale to the influence of a magnetic field. Do not weigh magnetic substances.

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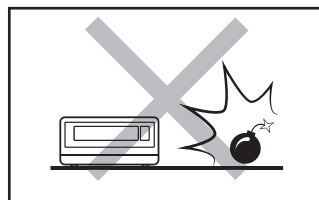Not intended for use within EX zones. Not intended for weighing explosive or flammable substances.

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
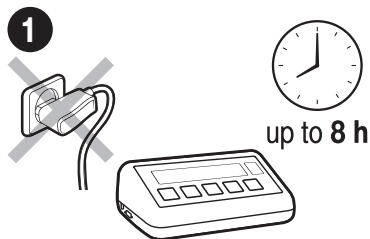
Prior to switching the device on let it reach room temperature (estimated stabilization time: 8 hours).

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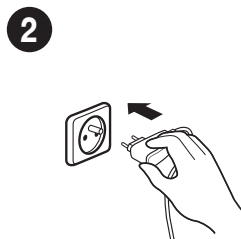
Connect the power supply to the mains.

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3.2. Indicator Start-Up

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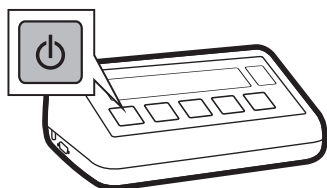Press ON/OFF key to start the indicator.

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The indicator is ready for operation!

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
4. PANEL AND SCREEN

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4.1. Keypad

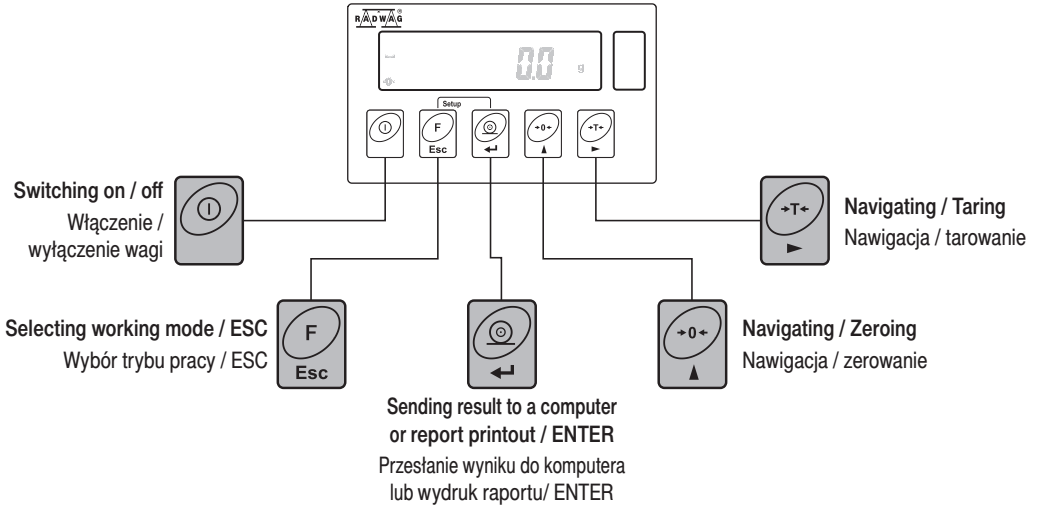Klawiatura

4.2. Home Screen

Ekran główny

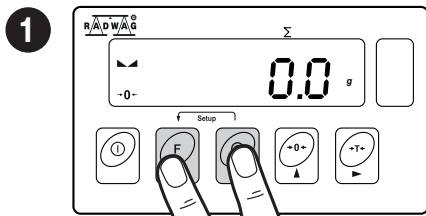
5. BASIC OPERATIONS

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
5.1. Working Mode Accessibility

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
Simultaneously press **Esc F** + **Enter** keys to enter the home screen.

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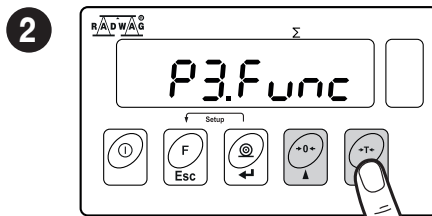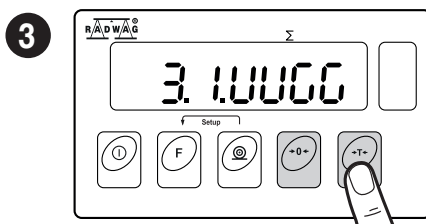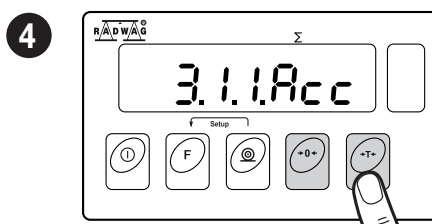
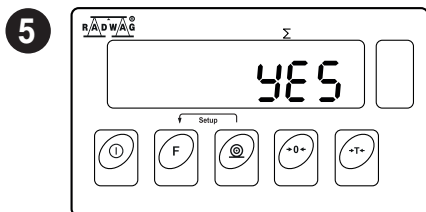
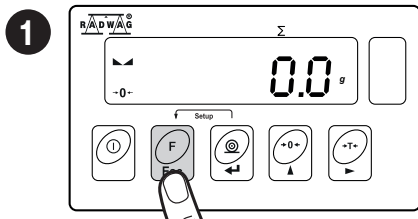
Follow the above procedure to configure other working modes.

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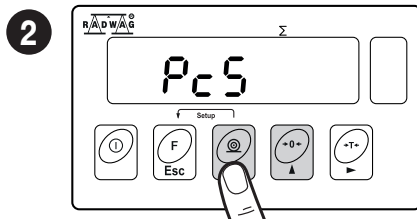
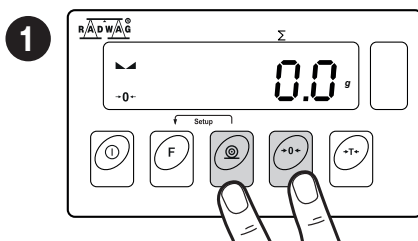
5.3. Unit Selection

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
When g (gram) is the main unit, you can select the following units: g (gram), ct (carat), lb (pound). *

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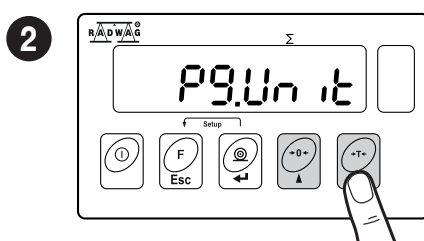
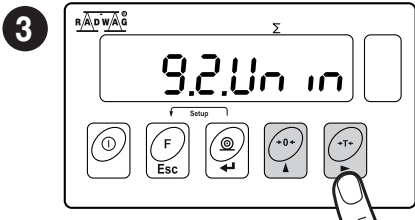
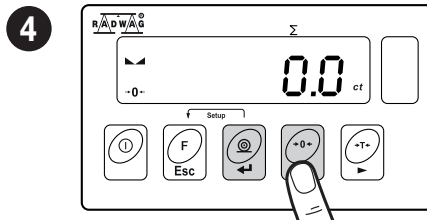
* Unit pound (lb) not available for verified scales.

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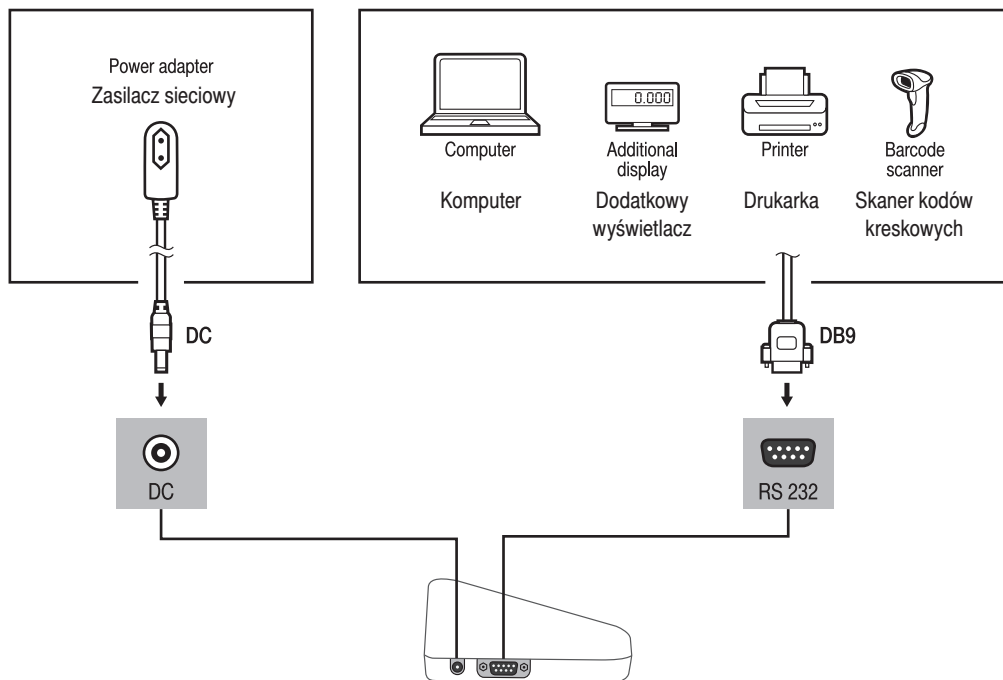
7. CLEANING AND MAINTENANCE

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
7.1. Notes and Warnings

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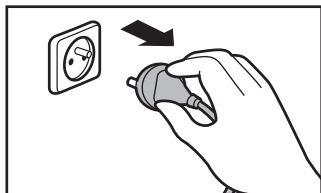
Prior to cleaning it is necessary to disconnect the scale from the mains. Unplug the connector from the electrical socket.

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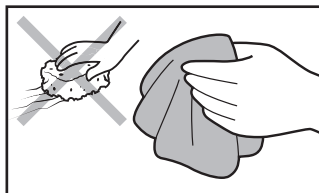For cleaning use only soft cloths made of microfiber. Do not use any abrasive cloths or any material that may scratch the surface!

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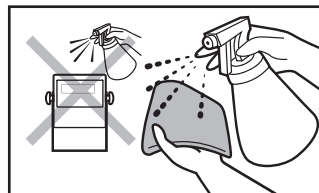Apply the cleanser onto the cloth first. Avoid applying the cleanser directly onto the device!

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Avoid using aggressive cleaning agents (e.g. solvents, cleansers containing chlorine, corrosive substances or bleach).

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Do not use cleansers containing abrasive substances or scouring preparations.

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Make sure that no dust or liquid enters the sockets and the inside of the indicator.

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